



**INSTRUCTIONS FOR SAMPLE COLLECTION AND SHIPMENT
FOR DUO ICS-KARYO TEST**








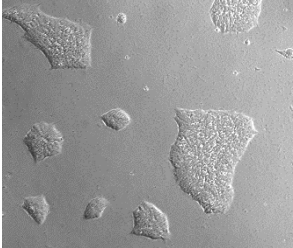
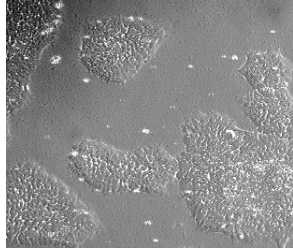

Please follow these instructions carefully to ensure proper sample reception and completion of the test.

Please do not send any sample to us without prior notice (1 week min.)

Samples should be received at Stem Genomics by Wednesday at the latest to guarantee timely processing of cells.

To be done at least one week before your shipment

Fill in the [Sample release form](#) carefully for all the samples to be tested and send it by e-mail in an Excel format to services@stemgenomics.com.

Method	Material	Shipment Temperature
<p>-Warm a cold pack to 37°C by placing it in an incubator few hours prior to the shipment.</p> <p>-Cultivate your cells for testing into two T25 flasks. The cell culture should reach about 40 and 60% confluency at the time of the shipment.</p>  <p>-Fill the T25 flasks to capacity with warmed media, tighten the cap and seal it with para-film. A 50 mL aliquot of media should be provided along with your cell cultures.</p>  <p>-Place samples and media into separate sealed bags.</p>  <p>-Place the warmed pack in the bottom of the package</p>  <p>-Pack the flasks and media securely to prevent them from shifting during transport.</p>  <p>-Place the printed version of the Sample release form completed inside the box.</p>	<p>For each sample (maximum 5 per shipment):</p> <p>One T25 flask at 40% confluency</p>  <p>+</p> <p>One T25 flask at 60% confluency</p>  <p>+</p> <p>50mL of media</p>	 <p>Room temperature (with the warmed pack)</p>



Shipment (the following instructions must be strictly followed)

Put the sample in the specific [UN 3373 packaging](#) "Biological substance, Category B" provided for free by most shipping companies.

Place the "This Side Up" labels on each side of the package.

Send the package [overnight](#) to:

Stem Genomics - Testing Service
IRMB - Hôpital Saint Eloi
80 avenue Augustin Fliche
34295 Montpellier - Cedex 5
France

Phone number and email address to give to the transporter
+33 985 036 160 - services@stemgenomics.com

Declared value: 10 € (or USD)
Description of Goods: samples
HS Code: 3822.00.0000 or 3822.00.00.90
The proforma invoice must contain the words "no commercial value" and "value for customs purposes only"

Contact services@stemgenomics.com with your shipping tracking number.