
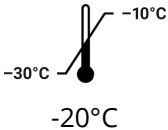





**INSTRUCTIONS FOR SAMPLE COLLECTION AND SHIPMENT
OF MYCO-DIGITAL TEST**

Sample	Method	Volume	Temperature	
			Storage	Shipment
 Cell culture supernatant	<ul style="list-style-type: none"> - The cells should be cultured without antibiotics for at least 3 days (this does not apply to Pen-step) - The culture medium needs to be in contact with the cells for minimum 24 h. - The cell culture should be at least 70% confluent. - Collect 500 µL of cell culture supernatant before passaging the cells and transfer into a safe-lock tube (DNase free) clearly labeled with your sample ID. - Boil supernatant at 95°C for 10 min 	<p>Tube 1: 500 µL of culture medium supernatant +</p> <p>Tube 2 (negative control): 500 µL of culture medium (without cell contact)</p>	 -30°C -10°C ≤72 h prior to shipment	 -20°C Room temperature


Fill in the [Sample release form](#) carefully for all the samples to be tested and [send it by e-mail in an Excel format to services@stemgenomics.com](#).



Shipment (The following instructions must be strictly followed)

The mention **UN 3373 "Biological substance, Category B"** is mandatory on the sample package, therefore your sample must be sent in four layers of packaging as follows:

- 1) Sample must be in a **leakproof tube**
- 2) Place the tube in a **first sealed bag with absorbent material** (for example paper towel). The maximum capacity allowed is 1L
- 3) Place the first sealed bag in a **second bigger and sealed bag**

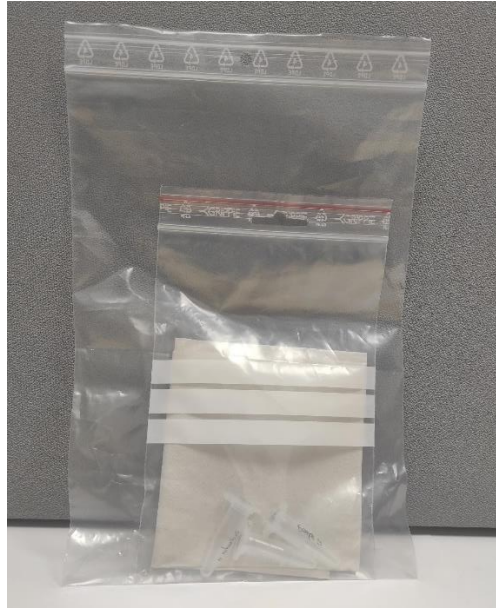

- 4) Place the sample in a **rigid packaging**, or a **bubble wrap envelop** or in the specific **UN 3373 packaging "Biological substance, Category B"** provided by most shipping companies free of charge.

Send it to:

Stem Genomics - Testing Service
IRMB - Hôpital Saint Eloi
80 avenue Augustin Fliche
34295 Montpellier - Cedex 5
France

Phone number and email address to give to the transporter:
+33 985 036 160 - services@stemgenomics.com

Declared value: 10 € (or USD)
Description of Goods: samples
HS Code: 3822.00.0000 or 3822.00.00.90
The proforma invoice must clearly indicate the mention "no commercial value" and "value for customs purposes only"