
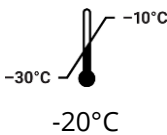








## INSTRUCTIONS FOR SAMPLE COLLECTION AND SHIPMENT OF MYCO-DIGITAL TEST


Sample	Method	Volume	Temperature	
			Storage	Shipment
 Cell culture supernatant	<ul style="list-style-type: none"> <li>- The cells should be cultured without antibiotics for <b>at least 3 days</b> (this does not apply to Pen-step)</li> <li>- The culture medium needs to be in contact with the cells for <b>minimum 24 h.</b></li> <li>- The cell culture should be at least <b>70% confluent.</b></li> <li>- Collect <b>500 µL</b> of cell culture supernatant before passaging the cells and transfer into a safe-lock tube (DNase free) clearly labeled with your sample ID.</li> <li>- Boil supernatant at <b>95°C for 10 min</b></li> </ul>	500 µL of culture medium supernatant	 <b>≤72 h prior to shipment</b>	 <b>Ambient temperature</b>
 Cells	<p>Dissociate cells (with enzymes, EDTA, etc.), centrifuge them (1200 rpm; 5 min), and discard the supernatant. Wash the cell pellet with PBS, centrifuge and discard supernatant.</p> <ul style="list-style-type: none"> <li>- Frozen cell pellets must be sent <b>on dry ice.</b></li> </ul>	Dry pellet <b>≥ 500,000 cells</b>		 <b>Dry ice</b>

Fill in the [Sample release form](#) carefully for all the samples to be tested and [send it by e-mail in an Excel format to services@stemgenomics.com](#).

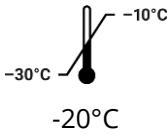**Shipment (The following instructions must be strictly followed)**

Your sample(s) must be sent in four layers of packaging as follows:

- 1) Sample(s) must be in a **leakproof tube**
- 2) Place the tube(s) in a **first sealed bag with absorbent material** (for example paper towel). The maximum capacity allowed is 1L.
- 3) Place the first sealed bag in a **second bigger and sealed bag**

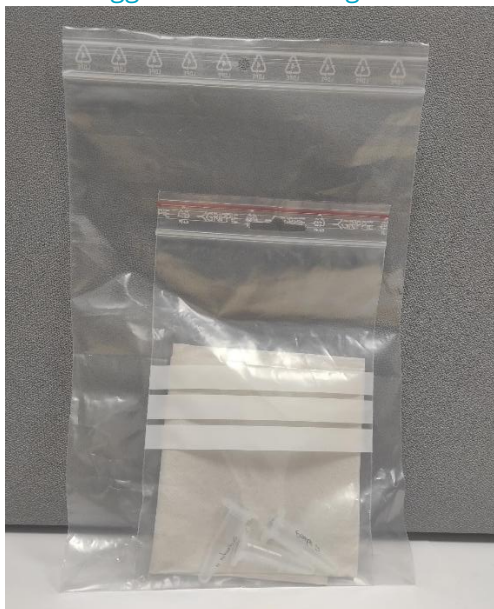

- 4) Place the sample in a **rigid packaging**, or a **bubble wrap envelop**.

Send it to:

Stem Genomics - Testing Service  
Cap Sigma  
1682 rue de la Valsière  
34790 Grabels  
France  
VAT n°: FR79841152077  
EORI n°: FR84115207700027

Phone number and email address to give to the transporter:

Phone : +33 420 900 201

Email : [services@stemgenomics.com](mailto:services@stemgenomics.com)

**Total declared value: 10 €**

**Description of Goods: samples, non-infectious, non-hazardous, not for medical purposes, for research purposes only. Goods packed in compliance with all applicable regulations. Not for human use.**

**HS Code: 3002.59 or 3002.5900 or 3822.00.0000 or 3822.00.00.90**

**The proforma invoice / commercial invoice must clearly indicate the mention "no commercial value" and "value for customs purposes only"**