
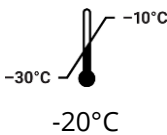








INSTRUCTIONS FOR SAMPLE COLLECTION AND SHIPMENT OF MYCO-DIGITAL TEST

Sample	Method	Volume	Temperature	
			Storage	Shipment
 Cell culture supernatant	<ul style="list-style-type: none"> - The cells should be cultured without antibiotics for at least 3 days (this does not apply to Pen-step) - The culture medium needs to be in contact with the cells for minimum 24 h. - The cell culture should be at least 70% confluent. - Collect 500 µL of cell culture supernatant before passaging the cells and transfer into a safe-lock tube (DNase free) clearly labeled with your sample ID. - Boil supernatant at 95°C for 10 min 	500 µL of culture medium supernatant	 -20°C ≤72 h prior to shipment	 Ambient temperature
 Cells	<p>Dissociate cells (with enzymes, EDTA, etc.), centrifuge them (1200 rpm; 5 min), and discard the supernatant. Wash the cell pellet with PBS, centrifuge and discard supernatant.</p> <ul style="list-style-type: none"> - Frozen cell pellets must be sent on dry ice. 	Dry pellet ≥ 500,000 cells	 -20°C	 Dry ice


Fill in the [Sample release form](#) carefully for all the samples to be tested and [send it by e-mail in an Excel format to services-us@stemgenomics.com](#).

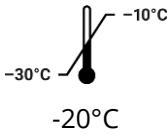

Shipment (The following instructions must be strictly followed)

Your sample(s) must be sent in four layers of packaging as follows:

- 1) Sample(s) must be in a [leakproof tube](#)
- 2) Place the tube(s) in a [first sealed bag with absorbent material](#) (for example paper towel). The maximum capacity allowed is 1L.
- 3) Place the first sealed bag in a [second bigger and sealed bag](#)

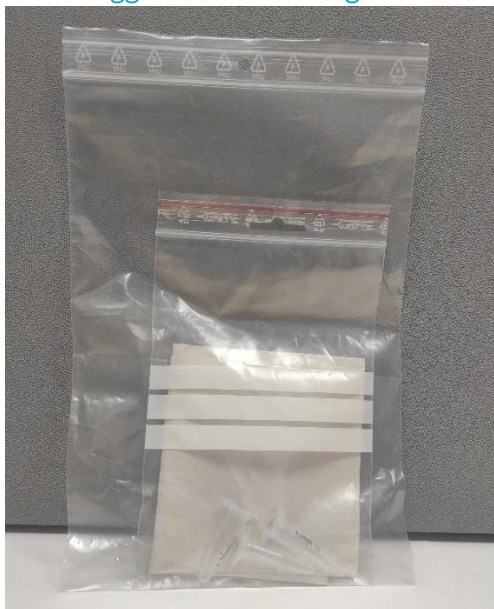

- 4) Place the sample in a [rigid packaging](#), or a [bubble wrap envelop](#).

Send it to:

Stem Genomics Inc. - Testing Service
400 Park Offices Drive, Room 104,
Research Triangle Park,
Durham, NC 27709
EIN n°: 38-4251349

Phone number and email address to share with
the transporter: services-us@stemgenomics.com
Phone: +1 (919) 638-5655

Total declared value: \$10

Description of Goods: samples, non-infectious, non-hazardous, not for medical purposes, for research purposes only. Goods packed in compliance with all applicable regulations. Not for human use.

HS Code: 3002.59 or 3002.5900 or 3822.00.0000 or 3822.00.00.90

The proforma invoice / commercial invoice must clearly indicate the mention "no commercial value" and "value for customs purposes only"